# MilkGuard Rapid test kit for Aflatoxin M1

## 1. Principle

This kit is based on the specific reaction of antibody-antigen and immunochromatography. Aflatoxin M1 in the sample competes for the antibody with the BSA-linked antigen coated on the membrane of the test strip. Then after a color reaction, the result can be observed.

## 2. Applications

This kit is used for rapid qualitative analysis of aflatoxin M1 in raw milk, Pasteurized milk or UHT milk.

## 3. Detection Limit (LOD): 0.5 µg/L(ppb)

#### 4. Kit components

- 8 wells and 8 test strips in one bottle, 12 bottles/ kit.
- microwell holder

- Kit insert
- Plastic pipette, 96pcs

#### 5. Operations

(1) Read the instructions carefully before experiment. Bring the test kit and samples to room temperature. Milk samples should be fully liquid without any agglomeration and deposition.

(2) Take bottles needed from the kit package, take out required wells and strips, and make proper marks. Please use these test strips within 1h. Seal the cap of the bottles.

(3) Take 200ul of the test samples into the wells, then repeatedly absorb for 5 times, mix the sample with the reagent in the wells completely. The mixture should be pink, and then start the timer.

(4) Incubate for 5min at room temperature(20~25°C), insert the test strips into the wells with the "MAX" end fully dipped in to the mixture.


(5) Incubate for 5min at room temperature(20~25°C) again. Take out the strip; judge the result according to Part 6.

#### 6. Result

**Negative(-)** : Line T and Line C are both red, and Line C is less stronger than Line T.

**Positive**(+) : Line C is red,Line C is stronger than Line T or Line C is the same color as Line T(aflatoxin M1>200ppt).

**Invalid**: Line C has no color, which indicates the strips are invalid. In this case, please

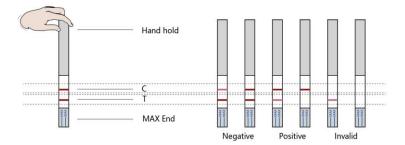

read the instructions again, and redo the assay with new strip.

Note: If the result of the strip needs to be recorded, please cut the foam cushion of the "MAX" end, and dry the strip, then keep it as file.

#### 7. Specificity

The kit shows negative with  $500\mu g/L$  of sulfamethazine, norfloxacin, lincomycin, spectinomycin, gentamicin, streptomycins. And shows positive with  $10\mu g/L$  of Aflatoxin G2,  $2\mu g/L$  of Aflatoxin B1, B2, G1.

#### 8. Storage

2-8°C in cool dark place, do not freeze. The kit will be valid in 12 months. The lot number and expired date are printed on the package.


#### 9. Notice for operations

(1) Please do the assay following the instruction, don not touch the membrane of the strip.

Beijing Kwinbon Biotechnology Co., Ltd

Add: No.8, High Ave 4, Huilongguan International Information Industry Base, Changping District, Beijing 102206, P. R. China Tel: +86-10-80700520/1 / 2 / 3. ext8811 Fax: +86-10-62711547 Web : <u>http://www.kwinbon.com</u> Email:xingyue@kwinbon.com P 1/2

- Please seal the bottle after taking out required strips. (2)
- This strip is used for only once; please do not use it repeatedly. (3)
- This kit is only for screening test, positive result should be further confirmed with other method. (4)

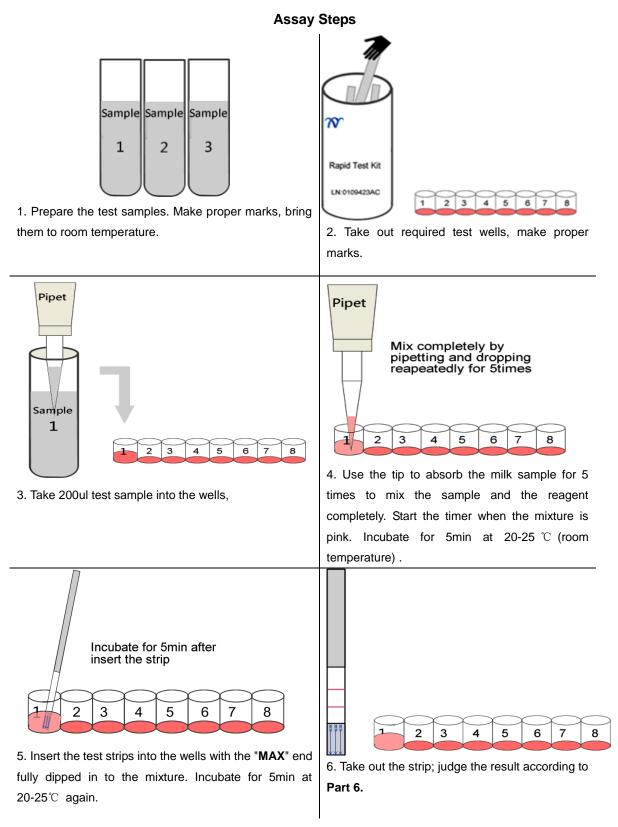

Beijing Kwinbon Biotechnology Co., Ltd Add: No.8, High Ave 4, Huilongguan International Information Industry Base, Changping District, Beijing 102206, P. R. China Tel: +86-10-80700520/1 / 2 / 3. ext8811 Fax: +86-10-62711547 P 2/2 Web : <u>http://www.kwinbon.com</u> Email:xingyue@kwinbon.com